



# Human TNF- $\alpha$ ELISA Test Kit Manual

Catalog #: 2105

*Reference #: 2105-01*

---

## TABLE OF CONTENTS

<b>GENERAL INFORMATION .....</b>	<b>2</b>
<i>Product Description .....</i>	<i>2</i>
<i>Procedure Overview.....</i>	<i>2</i>
<i>Kit Contents, Storage and Shelf Life .....</i>	<i>3</i>
<i>Required Materials Not Provided With the Kit .....</i>	<i>3</i>
<i>Sensitivity (Detection Limit).....</i>	<i>3</i>
<i>Specificity (Cross-Reactivity) .....</i>	<i>3</i>
<i>Warnings and Precautions .....</i>	<i>4</i>
<b>HUMAN TNF-<math>\alpha</math> ELISA TEST KIT PROTOCOL .....</b>	<b>5</b>
<i>Reagent Preparation.....</i>	<i>5</i>
<i>ELISA Testing Protocol.....</i>	<i>5</i>
<i>Human TNF-<math>\alpha</math> Concentration Calculations.....</i>	<i>7</i>
<b>TROUBLESHOOTING .....</b>	<b>8</b>
<i>No Color Development or No Signals with Standards.....</i>	<i>8</i>
<i>Low Optical Density (OD) Readings.....</i>	<i>8</i>
<i>High Background or High Optical Density (OD) Readings .....</i>	<i>8</i>
<i>High Intra-Plate or Inter-Plate Variance .....</i>	<i>9</i>
<i>One or More of the Standard Curve Points Are Out of Range.....</i>	<i>9</i>

*Human TNF- $\alpha$  ELISA Test Kit is intended for laboratory use only, NOT for clinical diagnostic use.*

## GENERAL INFORMATION

### Product Description

*Human TNF- $\alpha$  ELISA Test Kit* is designed for quantitative determination of the concentration of human tumor necrosis factor alpha (TNF- $\alpha$ , also known as cachectin) in serum, plasma, and cell culture supernatant.

Human TNF- $\alpha$  is a 17.5 kDa polypeptide of 157 amino acids. It is a potent lymphoid factor which exerts cytotoxic effects on a wide range of tumor cells. This cytokine is a primary regulator of inflammatory responses. TNF- $\alpha$  is produced by macrophages, neutrophils, activated T and B lymphocytes, endothelial cells, smooth muscle cells, astrocytes, natural killer cells, lymphokine-activated killer cells, and some transformed cells. TNF- $\alpha$ , and LT- $\alpha$  (lymphotoxin cellular receptors (CD 120a and CD120b), and produce similar effects. TNF- $\alpha$  plays a critical role in the destruction of certain tumors and in the body's resistance to infection by inducing fever and activating macrophages. Over-production of TNF has been associated with cytotoxic effects such as cachexia.

The specifications of the *Human TNF- $\alpha$  ELISA Test Kit* are as follows:

Format: 96-well strip plate

Assay range: 8-512 pg/mL

Sensitivity: 8 pg/mL

Total Assay Time: 3 hour and 45 minutes

Sample Size: 100  $\mu$ L/well

Sample types: Serum, Plasma and Cell Culture Supernatant

Wavelength for plate reading: 450 nm

### Procedure Overview

The method is a solid phase sandwich ELISA (Enzyme-Linked Immunosorbent Assay). It utilizes a monoclonal antibody (capture antibody) specific for human TNF- $\alpha$  coated on a 96-well plate. Standards and samples are added to the wells, and any human TNF- $\alpha$  present binds to the immobilized antibody. The wells are washed and biotinylated monoclonal anti-human TNF- $\alpha$  antibody (detection antibody) is added. After a second wash, avidin-horseradish peroxidase (avidin-HRP) is added, producing an antibody-antigen-antibody sandwich. The wells are again washed and a substrate solution is added, which produces a blue color in direct proportion to the amount of human TNF- $\alpha$  present in the initial sample. The stop buffer is then added to terminate the reaction. This results in a color change from blue to yellow. The wells are then read at 450 nm.

## **Kit Contents, Storage and Shelf Life**

*Human TNF- $\alpha$  ELISA Test Kit* has the capacity for 96 determinations or testing of 40 samples in duplicate (assuming 16 wells for standards and negative controls). Return any unused microwells to the foil bag and reseal them with the desiccant provided in the original package. Store the kit components as recommended in the table below. The shelf life is 12 months when the kit is properly stored.

<b>Kit Contents</b>	<b>Amount</b>	<b>Storage</b>
Capture Ab-coated Microtiter Plate	1 x 96-well plate (8 wells x 12 strips)	2-8°C
Human TNF- $\alpha$ Standard	2 x 20 $\mu$ L	-20°C
250X Detection Antibody	55 $\mu$ L	-20°C
250X Avidin-HRP	60 $\mu$ L	-20°C
5X Assay Diluent**	15 mL	2-8°C
20X Wash Solution**	28 mL	2-8°C
Stop Buffer**	20 mL	2-8°C
TMB Substrate**	12 mL	2-8°C

\*\* Components with the same XpressBio part numbers are interchangeable among XpressBio kits as long as they are used before their expiration dates.

## **Required Materials Not Provided With the Kit**

- Microtiter plate reader (450 nm)
- 10, 20, 100 and 1000  $\mu$ L pipettes
- Multi-channel pipette: 50-300  $\mu$ L (*Optional*)
- Distilled deionized water

## **Sensitivity (Detection Limit)**

8 pg/mL

## **Specificity (Cross-Reactivity)**

<b>Analytes</b>	<b>Cross-Reactivity (%)</b>
Human TNF- $\alpha$ :	100
Human IL-2:	<0.01
Human IL-3:	<0.01
Human IL-4:	<0.01
Human IL-5:	<0.01
Human IL-6:	<0.01
Mouse IL-1:	<0.01
Mouse IL-2:	<0.01
Mouse IL-3:	<0.01
Mouse IL-4:	<0.01
Mouse IL-5:	<0.01
Mouse IL-6:	<0.01
Mouse IL-7:	<0.01
Mouse IL-9:	<0.01
Mouse IL-10:	<0.01
Mouse IL-11:	<0.01
Mouse IL-12:	<0.01
Rat IL-2:	<0.01
Rat IL-4:	<0.01

## **Warnings and Precautions**

XpressBio strongly recommends that you read the following warnings and precautions to ensure your full awareness of ELISA techniques and other details you should pay close attention to when running the assays. More information can also be found in the Troubleshooting section. Periodically, optimizations and revisions are made to the kit and manual. Therefore, it is important to follow the version of the protocol included with the kit.

- The standard contains human TNF- $\alpha$ . Handle with particular care.
- Briefly centrifuge all tubes to collect the liquid at the bottom of the tube
- Do not use the kit past the expiration date.
- Do not intermix reagents from different kits or different lots. ANTIBODIES AND PLATES ARE KIT- AND LOT-SPECIFIC. Make sure that the standards, detection antibody, avidin-HRP, and diluent are mixed in correct volumes.
- Try to maintain a laboratory temperature of (20 – 25°C / 68 – 77°F). Avoid running assays under or near air vents, as this may cause excessive cooling, heating and/or evaporation. Also, do not run assays in direct sunlight, as this may cause excessive heat and evaporation. Cold bench tops should be avoided by placing several layers of paper towel or some other insulation material under the assay plates during incubation.
- Make sure you are using only distilled deionized water since water quality is very important.
- Incubations of assay plates should be timed as precisely as possible. Be consistent when adding standards to the assay plate. Add your standards first and then your samples.
- Add standards to plate only in the order from low concentration to high concentration as this will minimize the risk of compromising the standard curve.
- Always refrigerate plates in sealed bags with a desiccant to maintain stability. Prevent condensation from forming on plates by allowing them to equilibrate to room temperature (20 – 25°C / 68 – 77°F) before opening (plates provided in packaging contain desiccant).
- Be sure samples are properly stored. In general, samples should be refrigerated at 2-4°C for no more than 1-2 days. Freeze samples to -20°C or colder if they need to be stored for a longer period. Frozen samples can be thawed at room temperature (20 – 25°C / 68 – 77°F) or in a refrigerator before use.
- Human TNF- $\alpha$  in plasma, sera, and cell culture supernatant can be measured directly using this kit without extraction.

*XpressBio makes no warranty of any kind, either expressed or implied, except that the materials from which its products are made are of standard quality. There is no warranty of merchantability of this product, or of the fitness of the product for any purpose. XpressBio shall not be liable for any damages, including special or consequential damage, or expense arising directly or indirectly from the use of this product.*

## HUMAN TNF- $\alpha$ ELISA TEST KIT PROTOCOL

### Reagent Preparation

**IMPORTANT:** All reagents should be brought up to room temperature before use (1 – 2 hours at 20 – 25°C / 68 – 77°F). Make sure you read the “Warnings and Precautions” section on page 3. Solutions should be prepared just prior to ELISA test. Briefly centrifuge all tubes to collect the liquid at the bottom of the tube. All reagents should be mixed by gently inverting or swirling prior to use. Prepare volumes that are needed for the number of wells being run. Do not return the reagents to the original stock tubes/bottles. Disposable reservoirs are recommended to minimize the risk of contamination.

#### 1. Preparation of 1X Assay Diluent

Mix 1 volume of 5X Assay Diluent with 4 volumes of distilled water. In general, 50 mL of 1X Assay Diluent is sufficient for one whole plate.

#### 2. Preparation of Standards

Add 10  $\mu$ L of Standard solution to 1 mL of 1X Assay Diluent to prepare 512 pg/mL. Make serial 2X dilutions in the same diluent to make standards of 256 pg/mL, 128 pg/mL, 64 pg/mL, 32 pg/mL, 16 pg/mL., and 8 pg/mL. Use 1X Assay Diluent as negative control. Discard the remaining Standard solution after use.

#### 3. Preparation of 1X Wash Solution

Mix 1 volume of the 20X Wash Solution with 19 volumes of distilled water.

#### 4. Preparation of 1X Detection Antibody

Mix 1 volume of 250X Detection Antibody with 249 volumes of 1X Assay Diluent.

#### 5. Preparation of 1X Avidin-HRP

Mix 1 volume of 250X Avidin-HRP with 249 volumes of 1X Assay Diluent.

### ELISA Testing Protocol

Label the individual strips that will be used and prepare working solutions of reagents as shown in the following example. Adjust the total amount as needed for number of wells that will be used.

**Note:** Briefly centrifuge the standard, detection antibody & HRP tubes before opening to collect the liquid at the bottom of the tube

Component	Volume per Well	24 Wells
1X Detection Antibody	100 $\mu$ L	2.4 mL
1X Avidin-HRP	100 $\mu$ L	2.4 mL
1X Wash Solution	3.0 mL	72 mL
Stop Buffer	100 $\mu$ L	2.4 mL
TMB Substrate	100 $\mu$ L	2.4 mL

1. Add 100  $\mu$ L of 1X Assay Diluent to duplicate wells to serve as negative controls. Add 100  $\mu$ L of each dilution of Human TNF- $\alpha$  Standards in duplicate into different wells. (*⚠ Add standards to plate only in the order from low concentration to high concentration*).
2. Add 100  $\mu$ L of each sample in duplicate into different sample wells.
3. Incubate for 2 hours at room temperature (20 – 25°C / 68 – 77°F) ( *⚠ Avoid direct*

*sunlight and cold bench tops during the incubation. Covering the plate with aluminum foil while incubating is recommended).*


4. Aspirate all fluid from each well after the incubation and wash wells 3 times with 250  $\mu$ L of 1X Wash Solution per wash. It is not necessary to agitate the plate during the wash steps. Allow Wash Solution to remain in wells for 1-2 min during each wash step. After last wash, invert the plate and gently tap the plate on paper towels to remove residual fluid. ( $\emptyset$  *Perform the next step immediately after plate washings. Do not allow the plate to air dry between working steps).*
5. Add 100  $\mu$ L of 1X Detection Antibody into each well and incubate the plate for 1 hour at room temperature.
6. Aspirate and wash the plate 3 times with 250  $\mu$ L of 1X Wash Solution per wash. After the last wash, invert the plate and gently tap the plate on paper towels ( $\emptyset$  *Perform the next step immediately after the third wash. Do not allow the plate to air dry between steps).*
7. Add 100  $\mu$ L of 1X Avidin-HRP into each well. Incubate the plate for 30 minutes at room temperature (20 – 25  $^{\circ}$ C / 68 – 77  $^{\circ}$ F).
8. Wash the plate 3 times with 250  $\mu$ L of 1X Wash Solution per wash. After the last wash, invert the plate and gently tap the plate dry on paper towels ( $\emptyset$  *Perform the next step immediately after the third wash. Do not allow the plate to air dry between steps).*
9. Add 100  $\mu$ L of TMB Substrate and incubate for 15 minutes at room temperature (20 – 25  $^{\circ}$ C / 68 – 77  $^{\circ}$ F). Start timing the reaction immediately after adding the substrate to the last well. Mix the solution by gently rocking the plate manually during the first minute of incubation. ( $\odot$   $\emptyset$  *Do not put any substrate back into the original container which could lead to contamination. Any substrate solution exhibiting coloration is indicative of deterioration and should be discarded. Covering the plate with aluminum foil while incubating is recommended).*
10. Add 100  $\mu$ L of Stop Buffer in the order of adding TMB substrate to stop the enzyme reaction.
11. Read the plate as soon as possible following the addition of Stop Buffer. Read plate on a plate reader at 450 nm wavelength ( $\emptyset$  *Before reading, wipe the bottom of the plate with a lint-free tissue to remove any moisture or fingerprints that could interfere with the reading.*)

## **Human TNF- $\alpha$ Concentration Calculations**

A standard curve can be constructed by plotting the adjusted average absorbance for each reference standard against its concentration in pg/mL on a logarithmic scale as shown in the figure below. The adjusted average absorbance is obtained by subtracting the average of the negative control absorbance from the average of each observed absorbance.

To determine the corresponding concentration of human TNF- $\alpha$  in pg/mL in samples, plot the adjusted absorbance value for each sample on the standard curve. The figure to the right shows a typical human TNF- $\alpha$  standard curve.

A special program with Excel functionality, *Cytokine ELISA Analysis Program in Excel*, is available upon request to evaluate the test results.

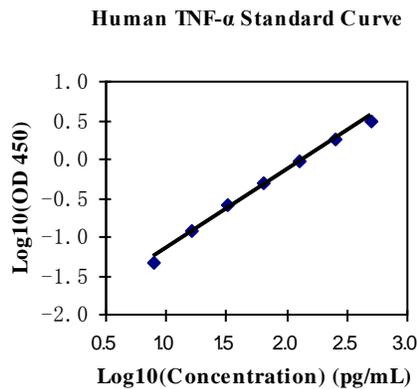

## TROUBLESHOOTING

### *No Color Development or No Signals with Standards*

Possible Causes	Recommended Action
<i>Reagents were used in the wrong order or a step was skipped.</i>	Follow the protocol carefully and repeat the assay.
<i>Wrong antibodies were used. Either Detection Antibody or Avidin-HRP was prepared incorrectly or has deteriorated.</i>	Make sure that the antibodies used are the ones that came with the kit. All antibodies are kit- and lot-specific. Make sure that the Detection Antibody, Avidin-HRP and diluent are mixed in correct volumes.
<i>TMB Substrate has deteriorated.</i>	Use a new set of TMB substrate. Note, if TMB substrate shows any color before use, it should not be used for the assay.

### *Low Optical Density (OD) Readings*

Possible Causes	Recommended Action
<i>Reagents were expired or mixed with a different lot number.</i>	Verify the expiration dates and lot numbers.
<i>Wash solution was prepared incorrectly.</i>	Verify that the solution was prepared as described in the protocol.
<i>Too many wash cycles were used.</i>	Make sure to use the number of washes per the protocol instruction.
<i>Incubation times were too short.</i>	Follow protocol and ensure accurate incubation time.
<i>Lab temperature was too low.</i>	Maintain the lab room temperature within (20 – 25°C / 68 – 77°F). Do not run assays under air conditioning vents or near cold windows.
<i>Reagents and plates were too cold.</i>	Make sure that the plates and reagents are brought up to room temperature.
<i>Reader was at wrong wavelength, or reader was malfunctioning.</i>	Make sure that the wavelength is set to 450 nm and read the plate again. Verify reader calibration and lamp alignment.
<i>Excessive kit stress has occurred.</i>	Check records to see how many times the kit has cycled from the refrigerator. Check to see if the kit was left at extreme temperatures for too long.
<i>Assay plates were compromised.</i>	Always refrigerate plates in sealed bags with a desiccant to maintain stability. Prevent condensation from forming on plates by allowing them to equilibrate to room temperature (20 – 25°C / 68 – 77°F) while in the packaging.

### *High Background or High Optical Density (OD) Readings*

Possible Causes	Recommended Action
<i>Poor quality water was used in wash solution.</i>	If water quality is questionable, try substituting an alternate source of distilled deionized water to prepare the wash solution.
<i>Substrate solution has deteriorated.</i>	Make sure that the substrate is colorless prior to addition to the plate.
<i>There was insufficient washing or poor liquid handling technique.</i>	Use the number of washes per the protocol instruction. Make sure that 250 $\mu$ L of wash solution is dispensed per well per wash. If you use a multichannel pipette or robotic liquid handling system, verify its performance; have the system repaired if any ports drip, dispense or aspirate poorly.
<i>Reader was malfunctioning or not blanked properly. This is a high possibility if the OD readings were high and the color was light.</i>	Verify the reader's performance using a calibration plate and check the lamp alignment. Verify the blanking procedure, if applicable, and reblank.
<i>Lab temperature was too high.</i>	Maintain the room temperature within (20 – 25°C / 68 – 77°F). Avoid running assays near heat sources or in direct sunlight.
<i>Reagents were intermixed, contaminated or prepared incorrectly.</i>	Ensure that the correct reagents were used, that working solutions were prepared correctly and that contamination has not occurred.

***High Intra-Plate or Inter-Plate Variance***

<b>Possible Causes</b>	<b>Recommended Action</b>
<i>Inconsistent time was taken when adding standards, reagents or samples within and/or between plates.</i>	Make sure that all materials are set up and ready to use. Use a multichannel pipette to add reagents to multiple wells whenever possible. Do not interrupt procedure while adding standards, reagents and samples.
<i>Multichannel pipette was not functioning properly.</i>	Verify pipette calibration and check that tips are on tight. Be sure all channels of the pipette draw and dispense equal volumes.
<i>There was inconsistent washing or poor liquid handling technique.</i>	Use the number of washes per the protocol instruction. Make sure that 250 $\mu$ L of wash solution is dispensed per well per wash. If you use a multichannel pipette or robotic liquid handling system, verify its performance; have the system repaired if any ports drip, dispense or aspirate poorly.
<i>Inconsistent incubation times occurred from plate to plate.</i>	Time each plate separately to ensure consistent incubation times.
<i>Pipette was inaccurate.</i>	Check the pipette calibration. Verify that pipette tips are on tight before use and that all channels draw and dispense equal volumes.
<i>Kit plates, reagents, standards and samples were at different temperatures.</i>	Make sure to allow sufficient time for kit plates, reagents, standards and samples to come to room temperature (20 – 25°C / 68 – 77°F). Larger volumes will require longer equilibration time. If using a water bath to hasten equilibration, make sure that it is maintained at room temperature; do not use a warm water bath to warm reagents, samples and kit standards.
<i>Reagents used were intermixed from different kit lots, or the kits were of different expiration dates.</i>	Carefully label each user-prepared reagent to make sure that the reagents are not intermixed. Kits with different expiration dates might generate different range of OD readings, however, the relative absorbance values will typically be comparable. In general, a value of less than 1.0 reading for the highest standard may indicate deterioration of reagents.

***One or More of the Standard Curve Points Are Out of Range***

<b>Possible Causes</b>	<b>Recommended Action</b>
<i>Standards were added in wrong order or recorded in wrong position.</i>	Make sure that the standards are applied and recorded correctly.
<i>Standards were contaminated or intermixed with other standards.</i>	Prepare a new set of standards. Always add standards to plate in the order from low concentration to high concentration.
<i>There was inconsistent washing or poor liquid handling technique.</i>	Use the number of washes per the protocol instruction. Make sure that 250 $\mu$ L of wash solution is dispensed per well per wash. If you use a multichannel pipette or robotic liquid handling system, verify its performance; have the system repaired if any ports drip, dispense or aspirate poorly.
<i>Inconsistent time was taken to add standards and reagents to plate.</i>	Make sure all materials are set up and ready to use. Add standards to plate only in the order from low concentration to high concentration at undisturbed constant pace. Use a multichannel pipette to add reagents to multiple wells simultaneously to increase consistency.
<i>Multichannel pipette was not functioning properly.</i>	Verify pipette calibration and check that tips are on tight. Be sure all channels of the pipette draw and dispense equal volumes.



XpressBio  
 503 Gateway Dr W  
 Thurmont, MD 21788  
 Tel: 301.228.2444  
 Fax: (301) 560.6570

Made in USA  
 info@xpressbio.com  
 www.xpressbio.com